

DNA-transfection of Madin-Darby canine kidney cells (MDCK II cell line) using Biontex K2® Transfection System

Lucie Navrátilová, Jana Mandíková, Milan Lázníček, František Trejtnar

Department of Pharmacology and Toxicology, Faculty of Pharmacy in Hradec Králové,
Charles University in Prague, Heyrovského 1203, 500 05 Hradec Králové, Czech Republic

Material and methods

Cell culture

The Madin–Darby canine kidney cell line (MDCK II) was cultured in 75 cm² cell culture flasks in the Dulbecco's modified Eagle's medium – high glucose (DMEM) containing 10 % fetal bovine serum in a humidified atmosphere containing 5 % CO₂ at 37 °C. Confluent cells were split twice a week using 0.25% trypsin/EDTA solution.

The MDCK II cells were seeded at a density 2×10^5 cells/well in 24-well plates. The transfection was performed the following day when the cells had reached a confluency of 90 %.

Cell transfection

The medium in each well was replaced with the fresh DMEM (0.5 ml) right before the transfection (before adding the K2® Multiplier).

2 h before DNA transfection 10 µl K2® Multiplier was added to the each well of cells and gently mixed by slightly moving of the plate.

Adequate amount of DNA and the K2® Transfection Reagent was added to the Opti-MEM® I Reduced Serum Medium (Gibco® by Life Technologies™) and gently mixed by pipetting up and down. Then the both solution was combined (DNA solution was added to the K2® Transfection Reagent solution), gently mixed and the mixture was incubated at room temperature for 20 min.

After incubation, 60 µl of DNA-lipid complex was added to the each well and gently mixed by shuffling of the plate. The cells were returned back to the incubator (5% CO₂, 37 °C) and incubated for 12 h or 24 h, respectively.

Cell line	K2® Multiplier [µl]	DNA [µg]	DNA/ K2® Transfection Reagent	Opti-MEM® [µl] for DNA + K2® Transfection Reagent
MDCK II	10	0.8	1:4	30+30


The transfection medium was replaced with the fresh DMEM after 12 h of the incubation and the results were analyzed after 24 h after the transfection. And the transfection medium was replaced with the fresh DMEM after 24 h of the incubation and the results were in this case analyzed after 48 h after the transfection.


Evaluation of the transfection

The success rate of the transfection after 24 h and 48 h was observed and compared with the cells transfected by the Lipofectamine® 2000 Transfection Reagent (Invitrogen™) using 0.8 µg DNA/well in the ratio DNA:Lipofectamine® 2000 1:2, 1:2.5 and 1:3. As a control served the MDCK cells transiently transfected with the empty vector (using the Lipofectamine® 2000 Transfection Reagent).

The MDCK cells were transfected with the OATP2B1 transporter and the accumulation study (24 h and 48 h after the transfection) was performed with the standard substrate of OATP2B1: radioactively labeled [³H]estrone 3-sulfate (ammonium salt) (E3S) (American Radiolabeled Chemicals).

Results

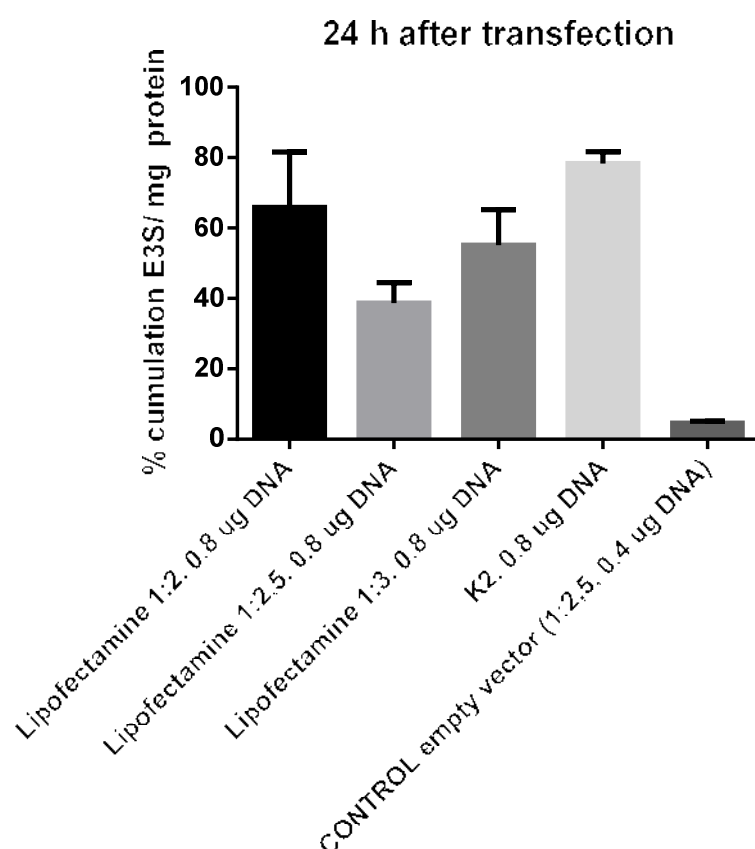

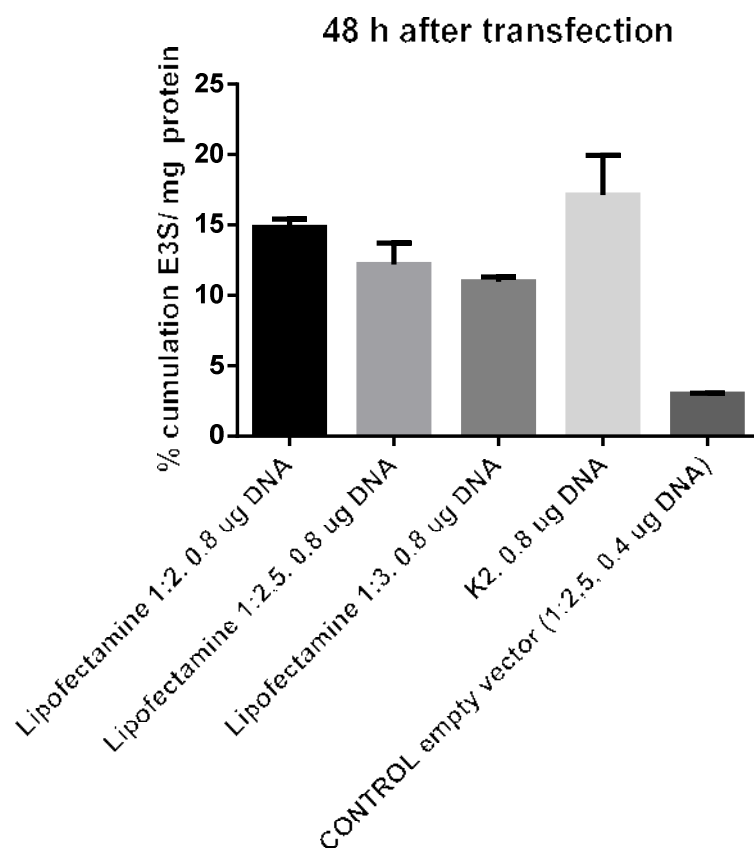

Conclusions

Transient transfection of MDCK II cells with the OATP2B1 expression construct was successfully performed using the K2® Transfection System. We observed **16.8** fold higher accumulation of [³H]estrone 3-sulfate in the cells transfected using K2® Transfection System versus the control after 24 h and **5.7** fold increase after 48 h. For the transient transfection of the MDCK II cells we obtained better results with the K2® Transfection System than the Lipofectamine® 2000 Transfection Reagent in the both cases. We achieved **1.2** fold increase of accumulation of [³H]estrone 3-sulfate using the K2® Transfection System versus using the most successful ratio of the Lipofectamine® 2000 Transfection Reagent (1:2, 0.8 µg DNA) after 24 h as well as after 48 h. These findings also confirm proper functionality of the MDCK II cell model transfected using the K2® Transfection System.