

Transient transfection of “DEP-1-HA in pNRTIS21” and “pNRTIS21” (negative control) into NIH3T3 using Lipofectamine (Invitrogen) and Metafectene Pro (Biontex)

Irene Weibrecht

(Prof. Ulf Landegrens group)

Uppsala University, Rudbeck Laboratory
Dept. Genetik och Patologi, Molecular Medicine
Dag Hammarskjölds Väg 20
75185 Uppsala, Sweden

Introduction:

Transient transfection of NIH3T3-cells with “hDEP-1-HA in pNRTIS 21” and “pNRTIS 21” (as negative control) should be done using Lipofectamine (Invitrogen) or Metafectene Pro (Biontex) for comparison. DEP-1-HA-expression and thereby transfection efficiency should then be investigated using immunofluorescence (IF). Dep-1-HA will be detected by mouse-anti-HA-antibody and staining with a donkey-anti-mouse-Cy3-labelled secondary antibody.

Materials and methods:

Transfection reagent: Lipofectamine (Invitrogen), Metafectene Pro (Biontex)

Cell-culture-medium: DMEM/F12 and DMEM (Sigma)

8-well-chamber slides (Lab-Tek II Chamber Slides)

Goat serum (Gibco)

Antibodies for IF: anti-HA-antibody (mouse-monoclonal, Nordic Biosite), donkey-anti-mouse-Cy3-labelled (Jackson Immunoresearch)

Experimental procedures / transfection protocol:

Lipofectamine:

-24 h prior to transfection 180,000 NIH3T3-cells were seeded per well of a 8-well chamber slide.

Day of transfection:

-Sol A: 0.15 µg DNA (DEP-1-HA in pNRTIS 21 and pNRTIS 21) were mixed with 30 µl DMEM/F12

-Sol B: 1 µl Lipofectamine (Lipofectamine: Invitrogen) was mixed with 30 µl DMEM/F12

-Sol A and Sol B were mixed and incubated at room temperature for 30 min (complex formation)

-The cells were washed with 300 µl/well DMEM/F12 and then 130 µl/well DMEM/F12 were added

-100 µl DMEM/F12 were added to the complexes and this solution was dropwise added to the cells

-The cells were incubated at 37°C, 5%CO₂ over night.

-The next morning the medium was changed to DMEM/F12 + 10% FCS + 50 U/ml penicillin/ 50µg/ml streptomycin and the cells were incubated for further 24 h to allow expression of the protein

Metafectene pro:

-16 h prior to transfection 320,000 NIH3T3-cells were seeded per well of a 8-well chamber slide.

Day of transfection:

-Sol A: 0.14 µg DNA (DEP-1-HA in pNRTIS 21 and pNRTIS 21) were mixed with 13.5 µl DMEM/F12 (mix by pipetting once)

-Sol B: 1.4 µl Metafectene Pro (Metafectene Pro: Biontex; Lot# AD2) was mixed with 13.5 µl DMEM/F12 (mix by pipetting once)

-Sol A was added to Sol B and incubated at room temperature for 20 min (complex formation)

-The medium of the cells was changed to 300 µl DMEM+10% FCS

-After 20 min the complexes were added dropwise to the cells.

-The cells were incubated at 37°C, 5%CO₂ for 48 h.

Fixation:

The slides were put on ice for 10 min, washed once with icecold 1x PBS and then fixed with 0.5% PFA in PBS on ice for 30 min.

The slides were washed with 1x PBS (icecold) and permeabilized using 70% EtOH (icecold) on ice for 30 min.

The slides were dried and stored at -20°C until use.

Immunofluorescence: The slides were blocked in 20% Goat serum, 0.05% Tween 20, 1x PBS at 37°C for 2 h. Primary mouse-anti-HA-antibody (100 ng/well) was applied in the same buffer and incubated at 4°C overnight. Staining with secondary donkey-anti-mouse-Cy3-antibody was done at room temperature for 1 h. Costaining of the nuclei was done using DAPI in the mounting medium. Pictures were taken with the 20x objective of an epifluorescence microscope.

Results and discussion:


See attached pictures. Transfection efficiency with lipofectamine was poor, the cell of which the picture was taken was among some ~5-10 cells/well which had been transfected. Transfection with Metafectene Pro was much better, here the picture is actually a representative picture for each field. In other words the transfection efficiency was much higher with Metafectene pro (~5%) compared to lipofectamine (~0.1%). The values are approximate.

Conclusion / summary:


For subsequent transfection of NIH3T3-cells Metafectene Pro will be used instead of Lipofectamine.

Appendix: Tables and/or figures:


METAFACTENE transfection with protein of interest
(DEP-1-HA)

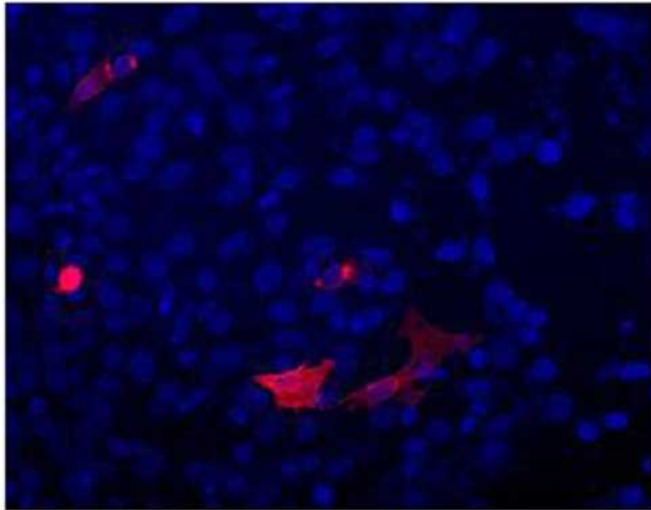

METAFACTENE transfection with empty vector
(negative control)

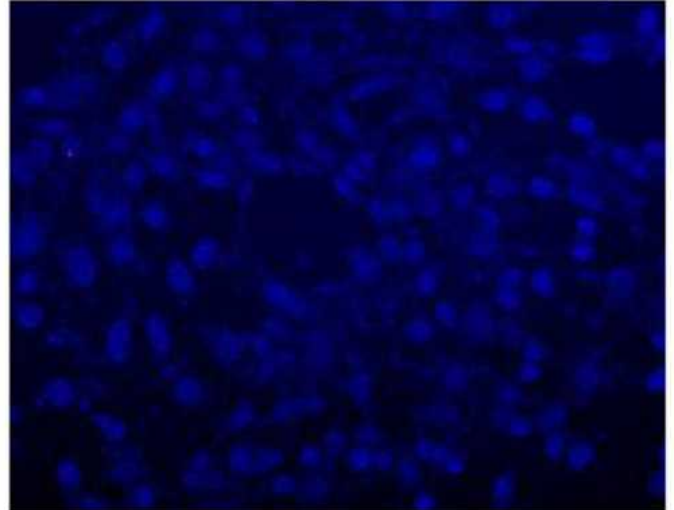

LIPOFECTAMINE transfection with protein of interest
(DEP-1-HA)

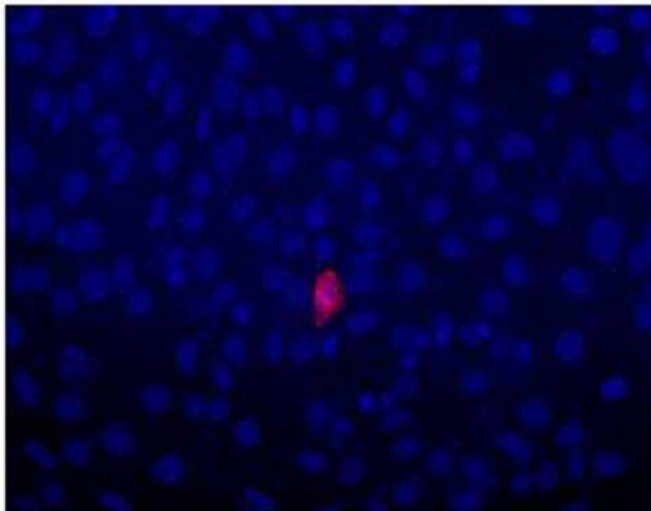

LIPOFECTAMINE transfection with empty vector
(negative control)

